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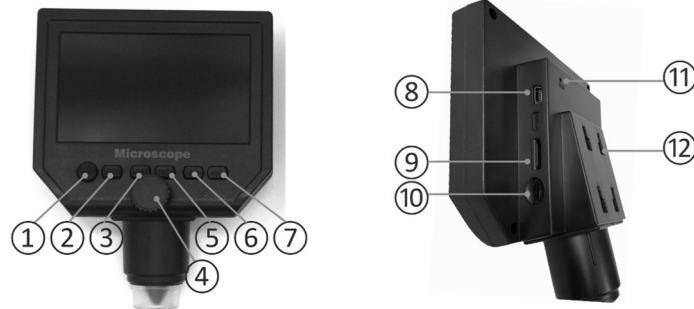
Warning: Please read this instruction manual carefully before using the instrument and keep it for future use.

A. SAFETY INSTRUCTIONS

1. The device is intended exclusively for private and not for commercial use. Only use this device as describe in this instruction manual.
2. Do not handle the unit or plug with wet hands.
3. When disconnecting the cable, always pull at the plug; do not pull on the cable itself.
4. If the cord is damaged, it must be replaced by the manufacturer or its service agent or a similarly qualified person in order to avoid an electric shock.
5. Never try to repair it by yourself. It must be repaired by the manufacturer or its service agent or a similarly qualified person in order to avoid an electric shock.
6. Always place your device on a dry surface.
7. Do not use outdoors.
8. Keep it away from heat sources such as radiators, to avoid deformation of plastic parts.
9. Do not use coarse or abrasive sponges/cloths or steel wool to clean the appliance.
10. This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capacities, or by persons lacking in experience or knowledge, unless they are supervised or instructed on the use of this apparatus by a person responsible for their safety. Children should be supervised to ensure they do not play with this appliance.

B. HOW TO USE IT

1) Charging the battery

Before using the microscope for the first time, connect the charger to the microscope and then plug it into an electrical outlet to charge the battery.

The symbol  will light up during the charging process and then change to the symbol  when the charge is complete (about 4 hours). Unplug the charger when the charge is complete to extend the life of the battery.

You will be able to use your device for about 6 hours before the next recharge. To prolong the life of the battery, remember to recharge it before the battery is completely discharged.

2) Inserting the SD card

Make sure the device is turned off

M 0600 / G600

Take a micro SD card.

Insert the micro SD card into the slot ⑨.

Without a memory card, the device cannot take pictures and videos directly.

Note: The device does not have a built-in memory. A memory card must be inserted into the camera before use.

3) Power on

- Press the Power button ① to turn on the device.

The device will automatically turn on in video mode, the logo  will be displayed on the top left of the screen.

- Press MENU button ② once, you can adjust the following settings by using the arrows ⑤ and ⑥ and OK button ⑦:

o Resolution: HD1280*720P or VGA 640x480

o Date: open / off

o Motion detection open / turn off (as soon as the camera detects motion it starts filming).

o Cyclic record: off / 3 minutes / 5 minutes / 10 minutes.

When the memory card is full, the first video is deleted and the new one is recorded automatically.

o Exposure value (EV): +2.0 (0) -2.0

- Press the MENU button ② a second time to adjust the following settings:

o Reticle: open / off

o Format: yes or no

o Language: choose one of the 16 languages

o Auto power off: off, 1 minute, 3 minutes,

o Screen savers: off, 1, 3, 5 minutes.

o Factory reset: Yes / No

o Light source frequency: 50HZ, 60HZ

o Time setting: year, month, day, hours, minutes, seconds.

o USB function: Memory / Webcam

- Press MODE button ③, the logo  will be displayed on the top left of the screen.

- Press MENU button ② once, you can adjust the following settings using the arrows ⑤ and ⑥ and OK button ⑦:

o Resolution: 5M, 3M, 2M, 1.3M, VGA

o Balance: automatic / sunlight / cloudy / tungsten / fluorescent

o Exposure: automatic / sunlight / cloudy / tungsten / Fluorescent / off 0 / -1/3 / -2/3 / -1 / -4/3 / -5/3 / -2

o Date: open / off

o Iso : automatic / 100 / 200

- When finished press MENU ② to exit the settings.

4) Operation

Attach the microscope to the flexible stand using the notches ⑫. Adjust the orientation of the objective to position the lens as close as possible to the object you wish to observe.

Place the object to be observed on the base and use the focus wheel ④ to get a sharp image.

Adjust the power of the LED light with the focus wheel ⑩.

In video mode press the OK button ⑦ to start video recording and press it again to stop it. If you have chosen motion detection mode, the camera will automatically record a video as soon as it detects motion.

In photo mode, press the OK ⑦ button to take a photo.

5) Viewing photos and videos

Press the MODE button ③ the logo  will appear on the top left of the screen.

Use the arrow keys ⑤ and ⑥ to select the file, and press OK ⑦ to play a video

C. TRANSFERRING RECORDINGS TO A COMPUTER

Turn off the device. Connect the device to the computer via the USB cable. Once the connection is established, a new dialog box appears on the computer. Photos and videos can be found in the "PHOTO" or "VIDEO" files in the mass storage and can be copied to the computer.

Note: Do not disconnect the cable or remove the memory card when copying photos/videos.

D. CARE AND MAINTENANCE

Clean the product only with a dry or slightly damp, lint-free cloth.

M 0600 / G600

Do not use abrasive cleaners to clean the unit.

E. FEATURES

- Materials: Plastic, 8 LEDs
- Power supply : Rechargeable lithium battery 2500 mAh, 3.7V, 9.25Wh (USB cable & AC adapter included)
- Magnification up to X 600 continuous - Resolution: 1080 / 720 / VGA
- Pixel HD 3.6 MP CCD
- Supports micro SD card up to 64 GB (card not included)
- Battery life : 6 h

The European directive 2012/19/EU about electronic and electric waste, requires that you can't throw away defective domestic appliance with common waste. Used device has to be collect separately to optimize the recuperation rate and the recycling of the materials to reduce the impact on the health and the environment.

Precision and precautions about the battery use

- Throwing away a used battery in the environment or with household waste pollutes and avoids the recuperation of reusable materials.
- It is important to limit the battery consumption and to respect the following rules:
 - Favor alkaline battery (last longer than saline battery) and when it's possible use rechargeable battery.
 - Leave used batteries and accumulators in adapted containers that you can find in some retailers. In that way, metals don't pollute the environment and the health.
- Battery has to be inserted by respecting the polarity show on the apparel and the battery. An incorrect positioning can damage the apparel, cause leaking and even provoke fire and the explosion of the battery.
- To assure a good functioning, battery has to be in a good condition. If there is a problem, put new battery.
- Never try to recharge battery not made for this. It could leak, warm up, cause fire or explode.
- Replace all the battery at the same time. Never mix up alkaline battery with saline battery.
- Used battery has to be removed.
- Also, remove the battery if you're not using the apparel for a long time.
- Never try to short-circuit the terminal of a battery.
- Never throw away the battery in fire because they can explode.
- Take out the accumulator from the apparel before reload it.
- We recommend that an adult watch if a child change the battery to respect those rules.
- If you swallow a battery, go immediately to your doctor or the hospital.
- For the best ongoing performance, we recommend using only fresh, high-quality, alkaline batteries.

Importé par / Importiert von / Geïmporteerd door / Imported by :
CAPTELEC, 59170 Croix - FRANCE

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